

Micro-Chamber/Thermal Extractor

Specification sheet

Covering: The four-chamber Micro-Chamber/Thermal Extractor (M-CTE250I).

The four-chamber Micro-Chamber/Thermal Extractor™ is a versatile, compact unit with four small cylindrical chambers that enables the sampling of chemicals released from products or materials at temperatures up to 250 °C.



1. System features

- **The system contains four inert-coated stainless steel microchambers** that allow surface or bulk emissions to be tested from up to four samples simultaneously.
- **A precision gas manifold** maintains a constant air/gas flow through each sample chamber.
- **No pump or mass flow controller is required** and near-ambient pressures are maintained in each test chamber.
- **Ambient or elevated temperatures** (up to 250 °C) can be used for sampling.
- **Total test time** (equilibration and vapour sampling), for all four samples, is normally between 15 and 30 min.
- **A conditioned sorbent-packed tube** is attached to each microchamber and a controlled flow of air/gas is passed through all chambers simultaneously.
- **Sampled sorbent tubes** are then loaded into a thermal desorption instrument (e.g. UNITY™) and analysed by TD-GC-MS.
- **A toggle valve option** is available that enables flows to individual chambers to be shut-off.

2. System specification

2.1 Dimensions and weight

- Height: 41.5 cm (16.3").
- Width: 16 cm (6.3").
- Depth: 52 cm (20.5").
- Weight: 15 kg (33 lb).

2.2 Microchamber dimensions

- Internal diameter: 6.4 cm.
- Depth: 3.6 cm.
- Available volume for bulk emissions testing: 114 cm³.
- Air volume above sample for surface emissions testing: 7.4 cm³.
- Exposed sample surface area for surface emissions testing: 24.63 cm².

2.3 Flow range (quoted for air or nitrogen)

- High: 50–500 mL/min through each chamber.
- Low: 10–70 mL/min through each chamber.

2.4 Gas specifications

- Gas types: clean air or nitrogen. Also compatible with helium, but flow rate range will vary.
- Gas pressure: 10–60 psig (0.69–4.15 bar).

2.5 Gas velocity at sample surface

- ~0.5 cm/s at 50 mL/min.
- ~5.0 cm/s at 500 mL/min.

2.6 Operational temperature range

- Ambient to 250 °C.
- Settable in 1 °C increments.

2.7 Ambient operating conditions

- Temperature: 15 °C to 30 °C.
- Relative humidity: 5 to 95% RH (non-condensing).

2.8 Power requirements

- 100–240 V, 50/60 Hz, 450 W maximum.

2.9 Safety and regulatory certifications

The instrument is designed and manufactured under a quality system registered to ISO 9001.

The instrument conforms to the following standards:

- International Electrochemical Commission (IEC):
 - 61010-1:2001.
 - 61010-2-010:2003.
 - 61010-2-081:2001.
- CAN/CSA C22.2 No. 61010-1 and UL 61010-1.

The instrument conforms to the following regulation on electromagnetic compatibility (EMC):

- IEC/EN 61326-1:2006.

For more information about our products and services, please visit www.markes.com.

Markes International shall not be liable for errors contained in this specification or for incidental or consequential damages in connection with the supply, performance or use of this equipment.

Markes International reserves the right to change information, descriptions and specifications within this document without notice.

Trademarks

Micro-Chamber/Thermal Extractor™ and UNITY™ are trademarks of Markes International.



Markes International Ltd

T: +44 (0)1443 230935 **F:** +44 (0)1443 231531 **E:** enquiries@markes.com

www.markes.com